

## NEULOG VOLTAGE LOGGER SENSOR GUIDE



### NeuLog voltage logger sensor NUL-201

The NeuLog voltage sensor can be used for any science experiment which utilizes voltage readings. It is used in the fields of Physics, Electronics, Chemistry, Biology, etc.

The sensor comes pre-calibrated so you can start experimentation right out of the box using this guide.

Just a few of the thousands of possible experimental subjects that can be studied with the NUL-201 sensor are: Van de Graaff generators, fuel cells, battery chemistry, human muscle physiology, parallel circuits, series circuits and electronics.

The voltage sensor's measurement units are:

Volts (V): The SI derived unit for electrical potential difference across two determined points.

### Quick start procedure:

#### PC or Mac Computer

#### Materials needed:

- NUL-201 Voltage Sensor
- USB-200 USB Module
- USB to mini USB cable (included with the USB-200 module)

Your voltage sensor needs to be connected to a USB-200 module. The USB-200 module then connects to a computer via a USB to mini-USB cable. Please note that you cannot plug the voltage sensor directly into the computer.

Resident PC software and browser based application can be downloaded for free at [www.neulog.com/Download/SoftwareandApplication.aspx](http://www.neulog.com/Download/SoftwareandApplication.aspx) as well as a full software user guide.

**Note:** Make sure not to download and install both types of software they will conflict on the computer.

#### Procedure:

1. Install the NeuLog software.
2. Connect the USB-200 module to the PC or Mac.
3. Connect the voltage sensor to the USB-200 module (they directly plug together). Please note no calibration is required for this sensor.
4. Open the NeuLog software.
5. A voltage sensor logo appears on the left side of the screen indicating that the probe has been automatically identified and you can begin experimentation.
6. If the voltage sensor is not automatically identified, click the "Search for sensors" icon to find the sensor.
7. Click the "On-line experiment" button; this will open a graph below.
8. Click on the "Module setup" button located on the voltage sensor icon in the module window to change the sensor's settings if need be.
9. Click on the "experiment setup" button to change the experiment settings if need be (experiment duration for example).
10. The voltage sensor will give a live reading in the box to the left of the screen while plugged in.
11. To run an experiment and collect data, click "Run experiment".
12. To end data collection early, click "Stop experiment".

#### Tablet, smart phone device

#### Materials needed:

- NUL-201 Voltage Sensor
- WiFi-201 WiFi Module
- BAT-200 Battery Module

Your voltage sensor needs to be connected to a WiFi-201 module. The WiFi-201 module will create a closed NeuLog WiFi network which will stream the NeuLog data to the device of your choosing. Once your device is wirelessly connected to the NeuLog network you can run experiments and collect data through a browser of your choosing.

#### Procedure:

1. Connect the voltage sensor directly to the left side of a WiFi-201 module (no wires required).
2. Connect a BAT-200 module to the right side of the WiFi-201 module.
3. Although not required, we recommend plugging the BAT-200 to an outlet using a USB to mini USB charger (such as a typical cell phone charger). The WiFi -201

## NEULOG VOLTAGE LOGGER SENSOR GUIDE



module will run for 15-75 minutes (depending on the sensor) without being plugged in.

- The WiFi-201 can be powered directly using a mini to USB cord and plugging it into your computer or a wall charger. Please note this function is only available on WiFi units with the USB icon on the cover.
- For further WiFi-201 instructions or the WiFi-201 quick start guide please visit: [www.neulog.com/Download/SoftwareandApplication.aspx](http://www.neulog.com/Download/SoftwareandApplication.aspx)

The WiFi-201 indicator lights will flash; take no action until the LED to the far left turns blue; this can take up to one minute.

- Take your tablet or smart phone, go to the Wi-Fi settings and select the NeuLog network which matches the WiFi Module ID found on the back of the WiFi-201 device. (NeuLogXXXX for example).
- Give your device about 20 seconds to connect to the WiFi-201.
- Once the device is connected, go to your browser and type the website [wifi201.com](http://wifi201.com) into the URL bar, then wait for 30-60 seconds.
- You will see a "Control mode" icon in the browser, click on this icon.
- The browser will then load a new screen and begin to auto detect the sensor(s); this can take a minute. (You can stop the search when the sensor is found.)
- If the browser does not auto detect the sensor(s), select "Search for sensors".
- Once the sensor is found, you will see an icon on the left side of the screen for the voltage sensor. The icon will display data in real time.
- To set the voltage sensor's settings, click on "Module setup" button located on the voltage sensor icon on the left side of your screen.
- To change the experimental settings, click the "Experiment setup" button (experiment duration for example).
- To run an experiment and collect data, select "On-Line" experiment, then select "Run".
- If you have a newer generation WiFi-201 (with the USB icon on the cover) this can be used as a USB-200, meaning you can tether the probes directly to the computer using the WiFi-201. To engage operation in this fashion, press 3 times on the panel key. Repeating this operation will return the unit to WiFi mode.

### Operation with NeuLog viewer

#### Materials needed:

- NUL-201 Voltage Sensor
- VIEW-101 Graphic Display Module
- BAT-200 Battery Module

#### Procedure:

- Connect the voltage sensor to the left side of the viewer.
- Connect the BAT-200 module to the right side of the viewer.
- The VIEW-101 will auto detect the sensor and you will see it appear on the left side of the screen.
- Once the sensor appears, it will be monitoring data in real time.
- To run an experiment and collect data, click the run button (little green person).

### Off-line experiments

(Off-line experiments are when you do not have a sensor connected directly to a computer, tablet, smartphone, or NeuLog viewer).

#### Materials needed:

- NUL-201 Voltage Sensor
- BAT-200 Battery Module

#### Materials needed to configure your offline experiment:

- USB-200 Module, WiFi-201 module or VIEW-101 Module
- USB to mini USB cable (included with the USB-200 module)

#### Procedure:

- Connect the voltage sensor directly to the left side of a charged BAT-200 module.
- When ready to collect data, press the "Start/Stop" button on the voltage sensor; a red light will turn on for the duration of the experiment.
- To change the experiment settings, first connect to your chosen device (PC, Mac, tablet, smart device) and select the "Module setup" button. (For more information on how to change the experiment settings, view the quick start procedure section for your chosen device in this document.)
- After the experiment has concluded (the "Start/Stop" button is pressed again or the data collection period ends) connect the sensor to your chosen device (you can find full instructions in each device's quick start procedure section).
- Open the NeuLog software.
- Click the "Off-line experiment" button.
- Click the "Load data from sensors" button.
- Select which experimental data to upload (5 experimental runs can be stored at one time on the NUL-201 voltage sensor).

## NEULOG VOLTAGE LOGGER SENSOR GUIDE



### Restoring sensor's factory default settings:

#### Procedure:

1. Connect the USB-200 module to a PC, Mac, tablet, or smart device.
2. Connect the voltage sensor to a USB-200 module (they directly plug together). Please note no calibration is required for this sensor.
3. Open the NeuLog software.
4. Click the "Tools" icon.
5. Click the "Restore sensor's factory defaults" icon.
6. If prompted to clear the graph, click either the "Off-line experiment" or "On-line experiment" button (whichever you are using).
7. Click "Clear experiment results".
8. Resume from step 4.

### Included with sensor:

- NeuLog NUL-201 Voltage Sensor Guide (this document).
- Red (positive probe) and black (negative probe) wires connected directly to the sensor's body.
- 4 mm red and black plugs for the positive and the negative wires.

Sensor's specifications	
Range and operation modes	±20 V
ADC resolution	12 bit
Accuracy	1%
Resolution	0.01 V
Max sample rate (S/sec)	100 for long duration 3000 for short duration

### Sensor's features:

- Fully digital data.
- Rugged plastic ergonomic case.
- 4 mm plugs for easy connectivity
- Push button switch for Start/Stop experiments in off line mode.
- LED indicator of experiment status (blinks while collecting data).

- Pre-calibrated sensing equipment.
- Note:** NeuLog products are intended for educational use.

### Videos and experiment examples:

- Videos, literature and other probes can be found at [www.NeuLog.com](http://www.NeuLog.com)  
<http://www.neulog.com/Products/VoltageLoggersensor.aspx>
- **Ohm's Law** - <http://www.neulog.com/Portals/3/Users/Experiment P-10 Ohm's Law Ver 3.4.4.pdf>

### Technical background:

The philosophy behind NeuLog's plug and play technology is based on each sensor's ability to store its own data due to an internal flash memory chip and micro-controller in each plastic NeuLog body. This technology allows the sensor to collect and then store the digital data in the correct scientific units (°C, °F, Lux, %, ppm, for example).

The sensor is pre-calibrated at the factory. The built-in software in the logger can be upgraded for free at any time using the provided firmware update.

The sensor is based on very fast, high input impedance differential operational amplifier with protected inputs.

### Maintenance and storage:

- Never submerge the NeuLog plastic body in any liquid.
- Do not allow liquid into the voltage sensor's body.
- After use, gently wipe away any foreign material from the voltage sensor.
- Store in a box at room temperature out of direct sunlight.

## NEULOG VOLTAGE LOGGER SENSOR GUIDE



### Warranty:

We promise to deliver our sensor free of defects in materials and workmanship. The warranty is for a period of 3 years from the date of purchase and does not cover damage of the product caused by improper use, abuse, or incorrect storage. Sensors with a shelf life such as ion selective probes have a warranty of 1 year. Should you need to act upon the warranty, please contact your distributor. Your sensor will be repaired or replaced.

Thank you for using NeuLog!



Flexible, simple, fast, forward thinking.

W: [www.neulog.com](http://www.neulog.com)

E: [info@neulog.com](mailto:info@neulog.com)

A: 850 St Paul Street, Suite 15, Rochester, NY 14605

P: 1.866.553.8536

V05162013